



Tel: (039058) 9762 -0
Fax: (039058) 9762 -22
E-Mail: info@shp-steriltechnik.de
Internet: www.shp-steriltechnik.de

Dear customer,

thank you very much for choosing SHP Steriltechnik AG product.

Your device is configured as **LABOKLAV MV**

to be able to operate device you need following connections:

- ☒ power supply: 3N 400V ~ , 50 Hz, 16A, CEE socket
- ☒ tap water (cold) – flow pressure $\geq 0,5\text{bar}$ equipped with cut-off valve, $\frac{1}{2}$ " outside thread
- ☒ de-ionized water – flow pressure $\geq 1\text{ bar}$ equipped with cut-off valve, $\frac{1}{2}$ " outside thread
if there is no de-ionized water connection available device has to be filled manually by user
- ☐ compressed air: 4-6 bar, pressure regulated, oil free and dry, output min. 200l / minute,
 $\frac{1}{2}$ " outside thread
- ☒ drainage: DN 50 reduced to $\frac{1}{2}$ " outside thread
(high of drainage at Laboklav 55-100 max. 300mm and at Laboklav 135-195 max. 600mm)

All connections have to be equipped with $\frac{1}{2}$ " outside thread

Distance from device connection to media connection should not be longer than 1,50 m
Flexible metal hoses for connection include in delivery (10 bar overpressure resistant, 1,5m long).

Please note that device installation can be done only if media connections are prepared.